



Flow Cytometry for Studying Exosomes and Extracellular Vesicles

SelectBIO Webinar

March 7, 2024

Giacomo Vacca, Ph.D.

Founder & CEO, Kinetic River

Prof. Christopher Ward

U. of Kansas Medical Center

Today's Webinar

Introduction to the *Delaware* Flow NanoCytometer®

G. Vacca

- Features and capabilities
- Results to date
- Configurations

Toward Flow-Based Rapid Diagnostics of Kidney Disease


C. Ward

- Polycystic Kidney Disease (PKD)
- Current methods and unmet needs
- Urine extracellular vesicles (EVs) as PKD biomarkers
- Noninvasive PKD dia/prognostics with ultrasensitive flow cytometry

Q&A



Kinetic River

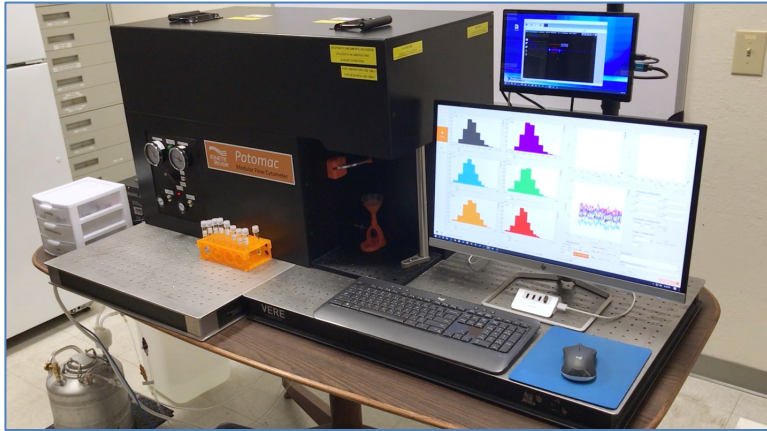

Mountain
View

California

- flow cytometers
- optical instrumentation
- design & development
- 24 issued patents, 7 pending
- past customers include: NIH-NCI, Stanford, UC Davis, Mt. Sinai, Millipore, Sony, NMSU, National Research Council (Italy)



Kinetic River Team



GIACOMO VACCA, PHD
Founder, President, & CEO



ALAN CHIN, PHD
Principal Scientist



ROSEMARY COATES, MBA
Advisor, Reshoring Institute



DAN ESTOQUE, BS
System Engineering
Consultant



RICH HANSON, MS
R&D Technician



KENNAN LANELUTTER, BS
Laboratory Technician



JKI
Software
Development Partner



ORIN LANEY, MS
Electronics Engineer
Consultant



HANNU LEHTIMÄKI, MS
Sr. Mechanical Engineering
Consultant



RICHARD MCKAY, PHD
Advisor, Full Spectrum
Scientific



SEAN MURPHY, MS
Advisor, SKMurphy




ASHLEY SLOAT, PHD
Advisor, Aurora Consulting



BONNIE SUN, BS
Office Administrator



CytoFlow Service
Distribution / Service
Partner, Europe

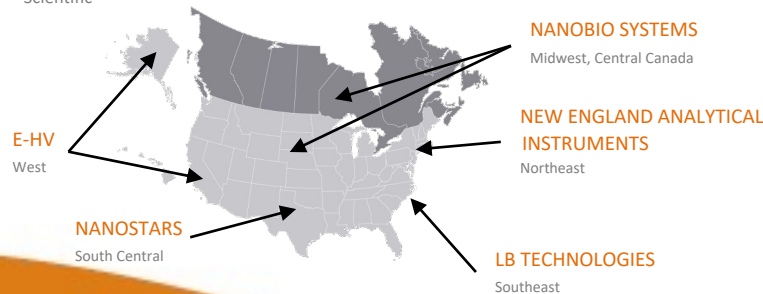


Delaware Flow NanoCytometer®



What Matters in EV Analysis?

- **Single-particle detection** (not bulk/averages)
- **Sensitivity** (smaller particles)
- **Multiparametric detection** (simultaneous biomarkers)
- **Throughput** (greater accuracy)
- **Flexibility** (measuring a wide range of particle sizes)
- **Stability** (confidence in results)
- **Ease of use** (focus on measuring, not tweaking)


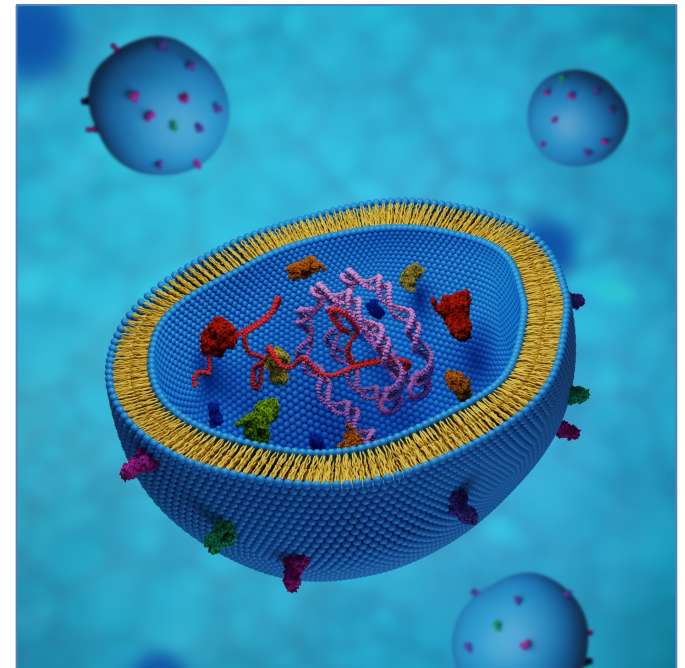


Delaware Flow NanoCytometer[®] for EV Analysis

- high resolution *and* sensitivity
- stable, pressure-driven flow
- up to 5 lasers
- up to 6 fluorescence channels
- measures EVs *and* cells



Delaware: Single-Particle Detection

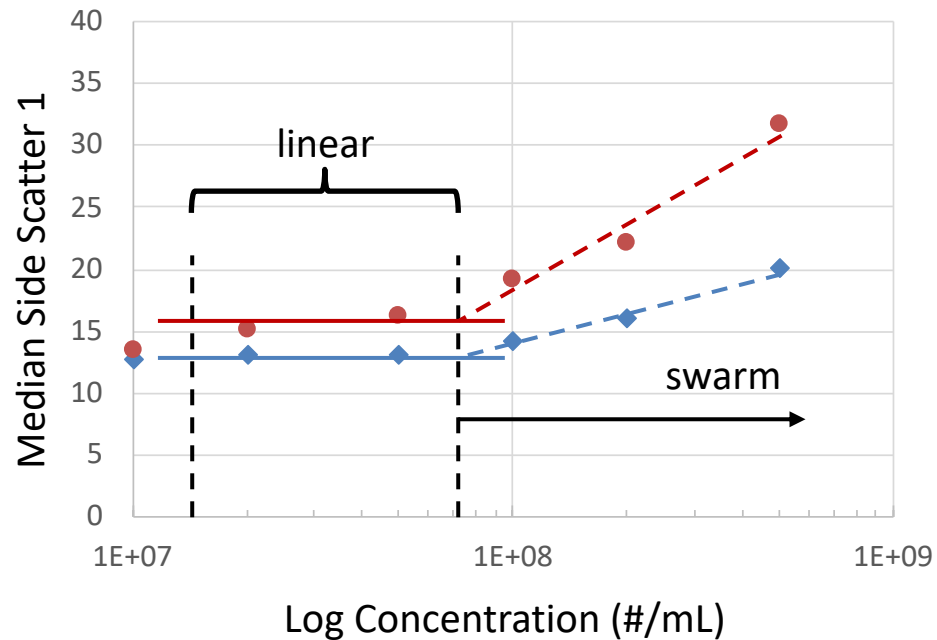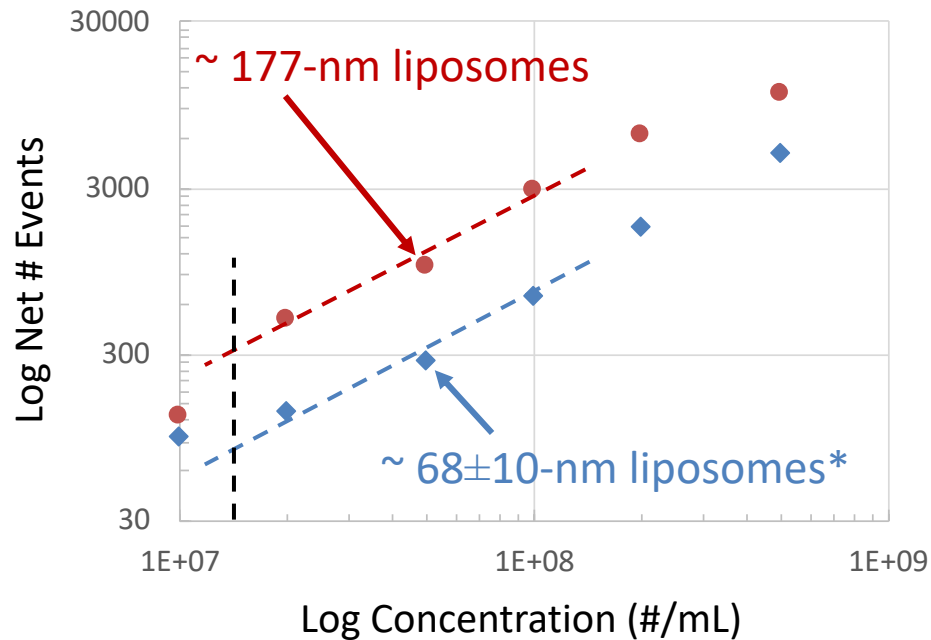


- liposome standards: Acoerela AcoLS

ACOERELA 

- sized by NanoFCM: 58 nm
- sized by ZetaView: 77 nm




Delaware: 60-nm Scattering Sensitivity (PS NPs)

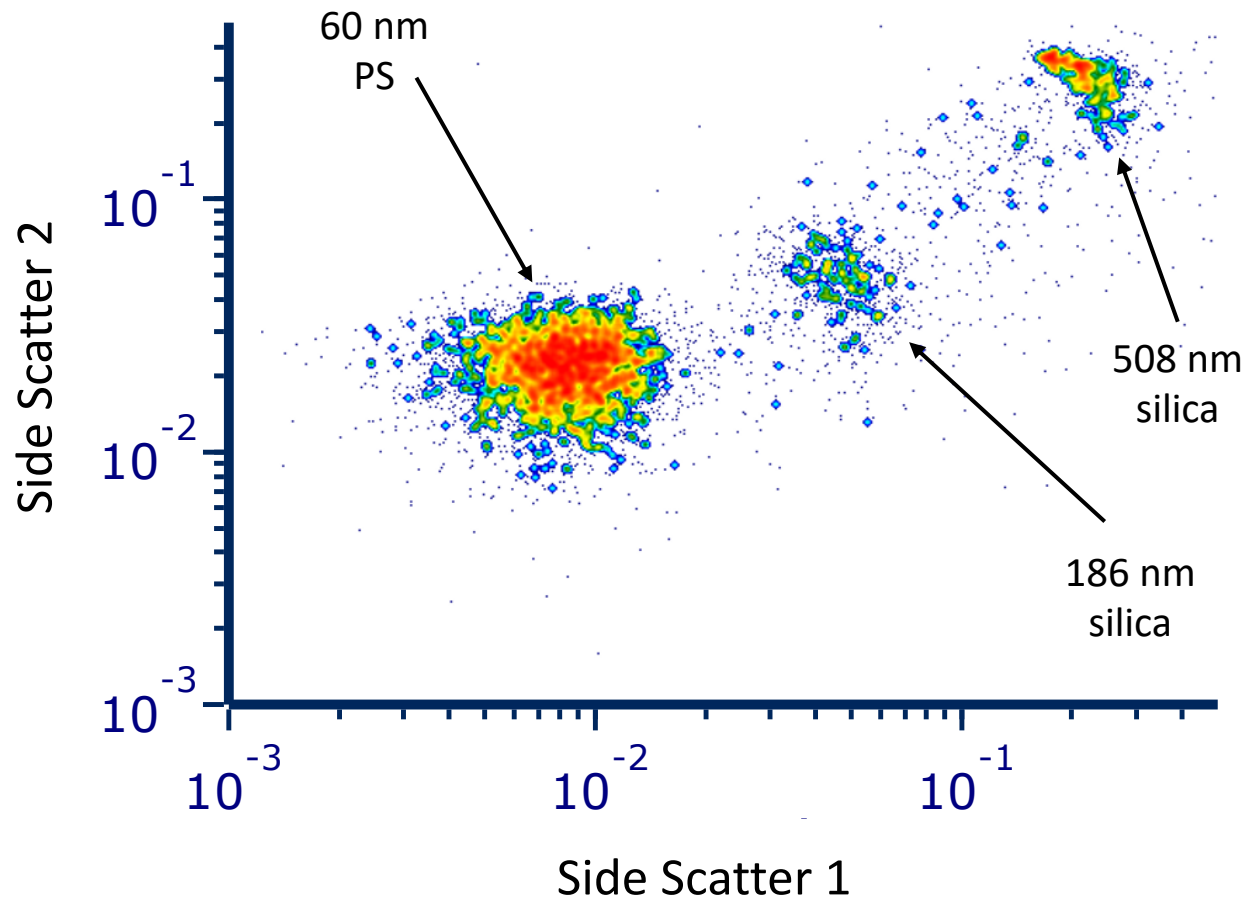


- silica: Alpha Nanotech
- polystyrene: Spherotech

Delaware: 28-nm Scattering Sensitivity (Au NPs)



102 nm Gold

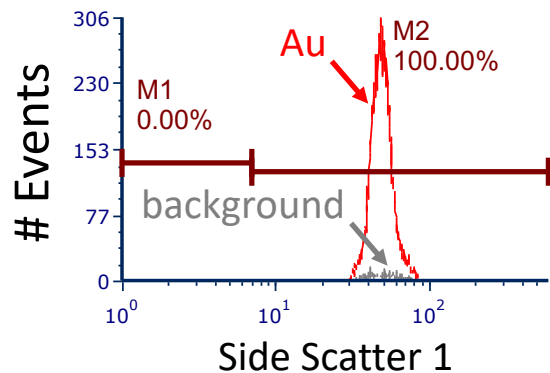

50 nm Gold

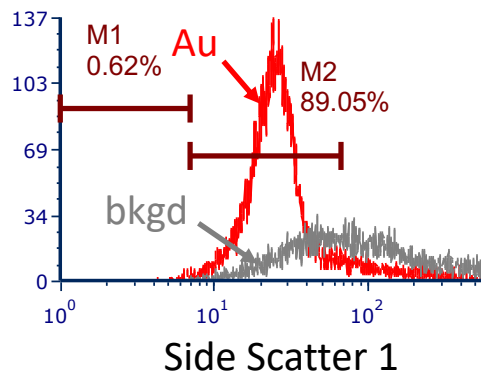


28 nm Gold

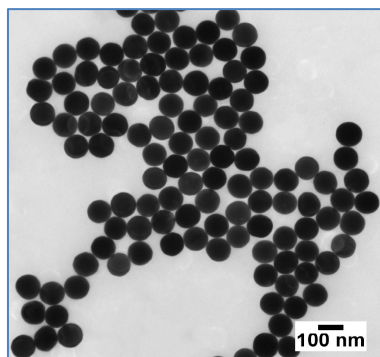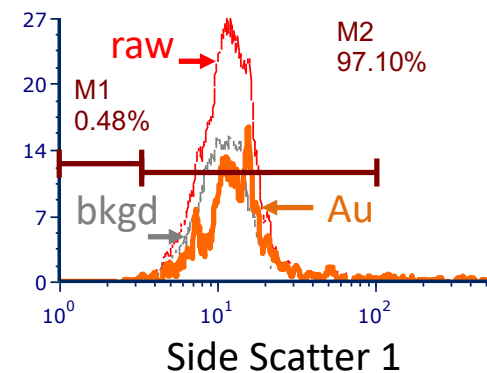


- Au NPs: nanoComposix
- sized by TEM

Delaware: 6-nm Resolution (2% CV)

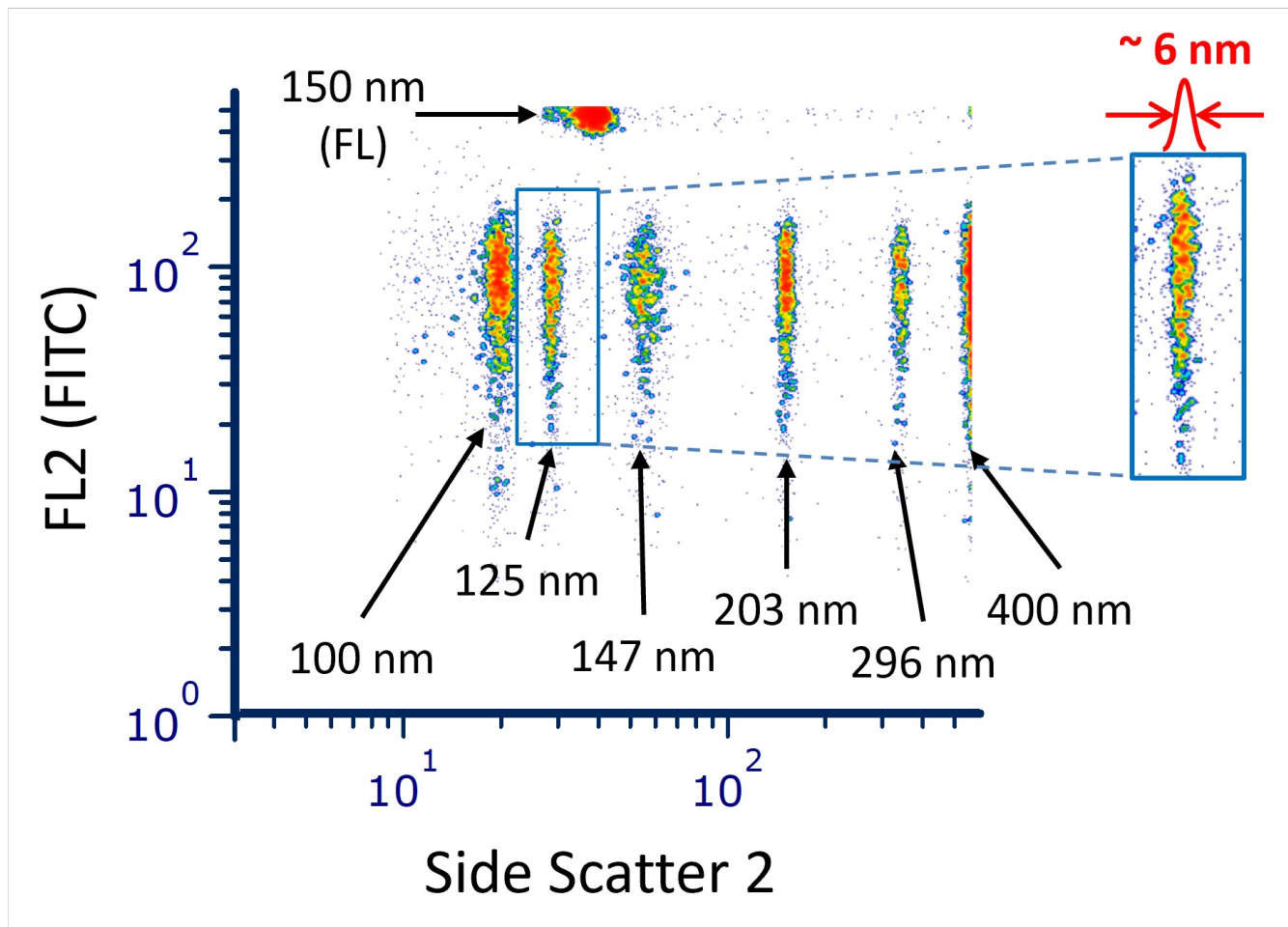

- nanoparticles:
Exometry Rosetta beads
- $CV = \sigma/\mu = FWHM / (2.36 * \mu)$


Delaware: Large Dynamic Range

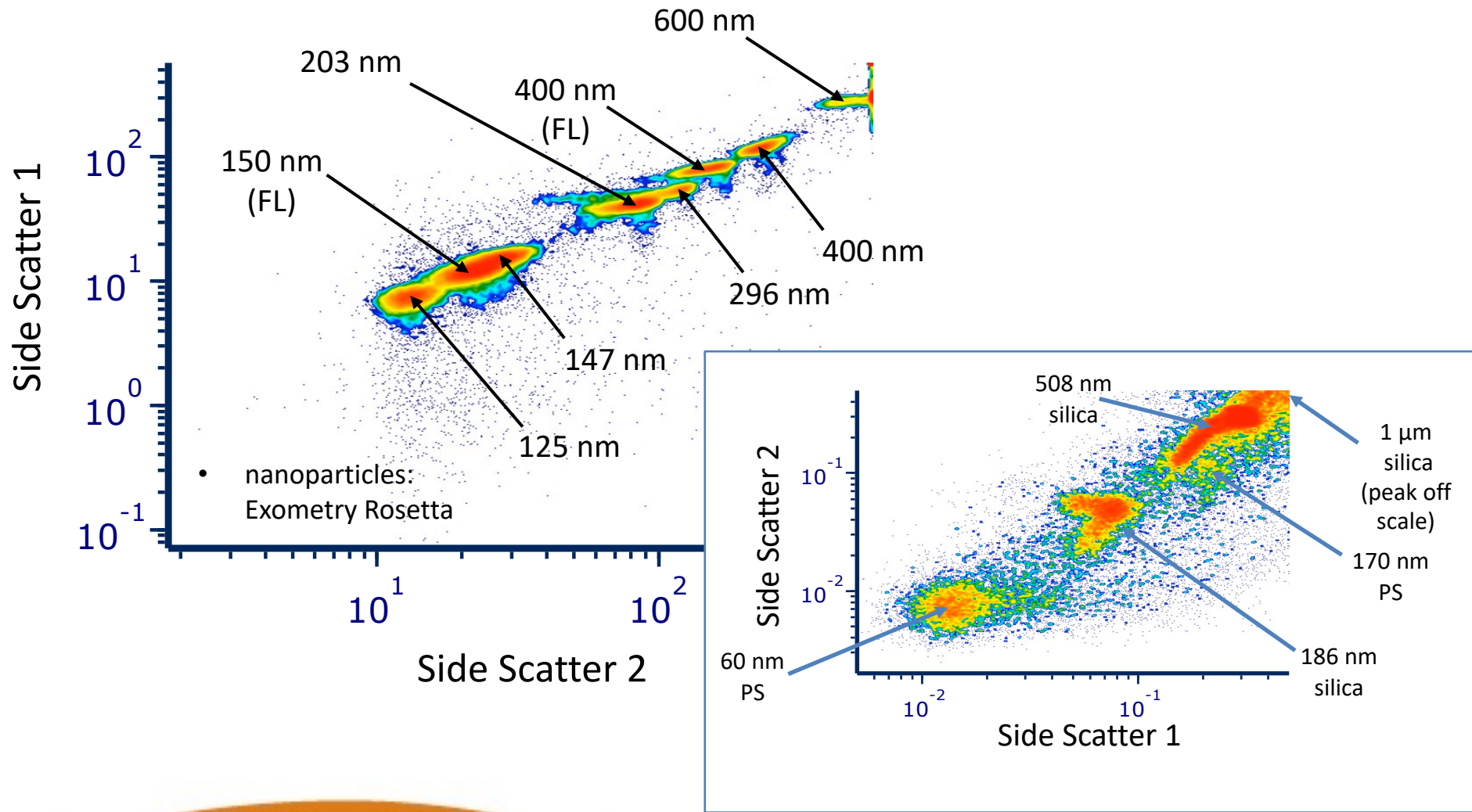




- silica: Alpha Nanotech
- polystyrene: Spherotech

Delaware: Human Cell-Derived EVs (Scattering)

HeLa EVs

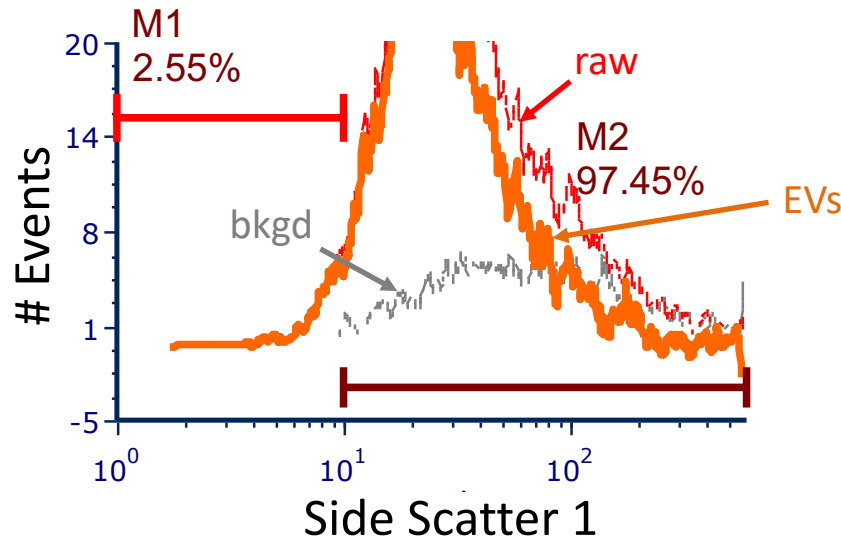


PC3 EVs

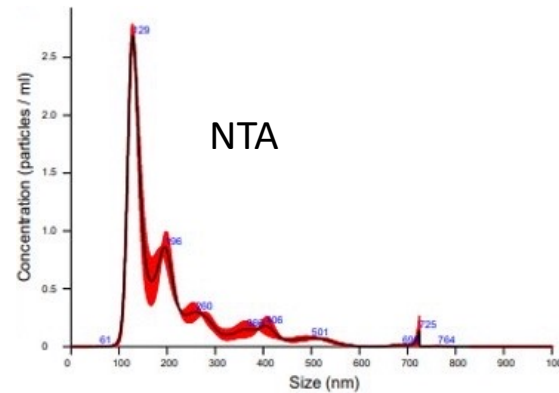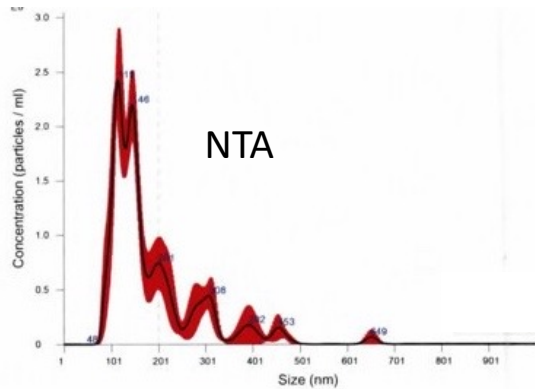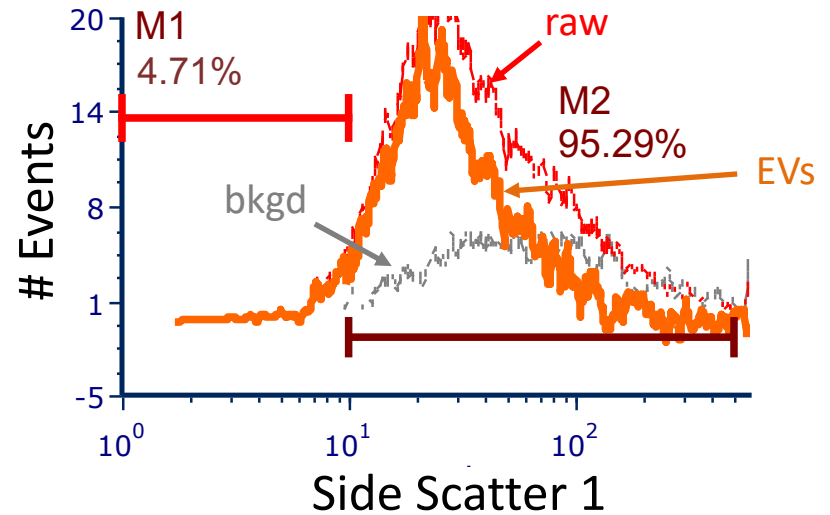








Delaware: Up to 6 Fluorescence Channels

- multiparametric phenotyping
- flow: complete, **simultaneous** colocalization, all the time

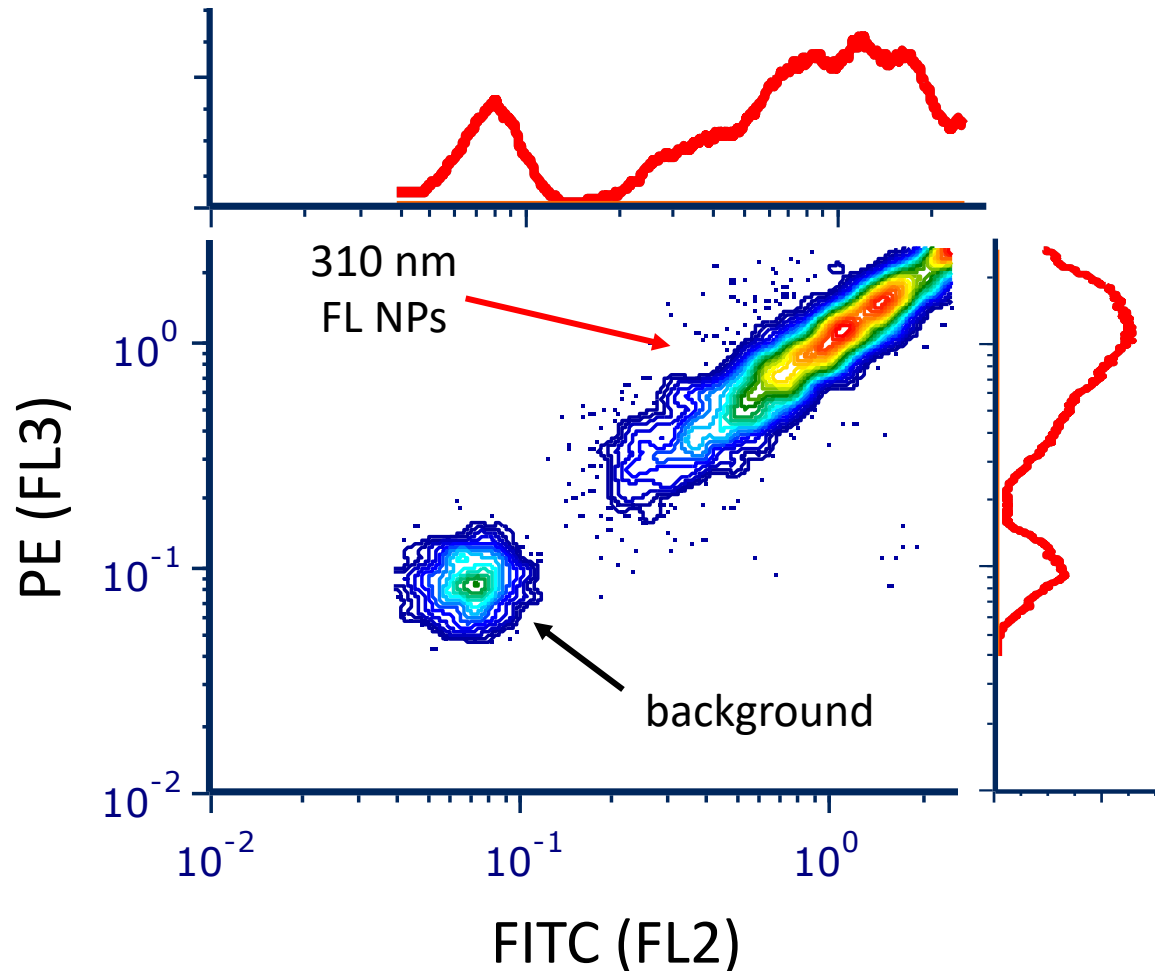

- fluorescent nanoparticles: Spherotech UltraRainbow 310 nm

Delaware: Human Cell-Derived EVs (Scattering, FL)

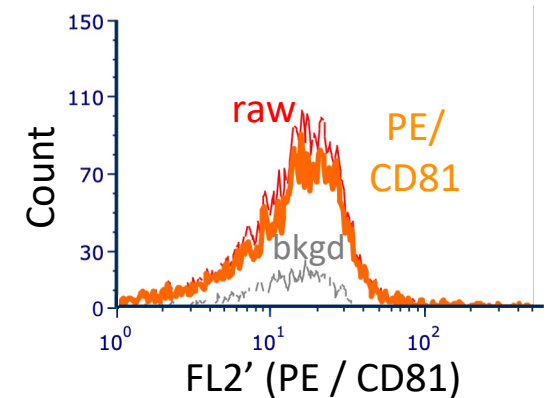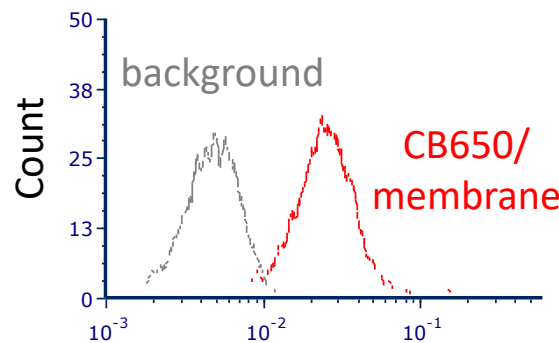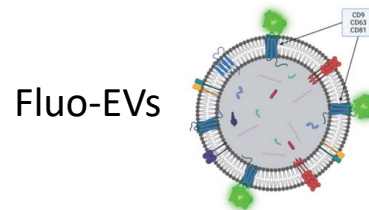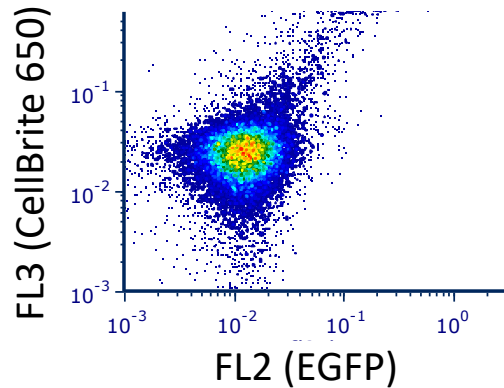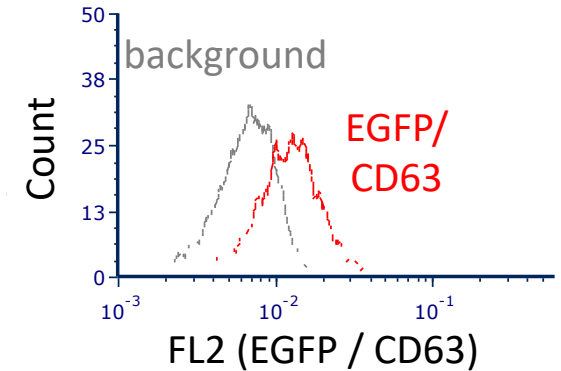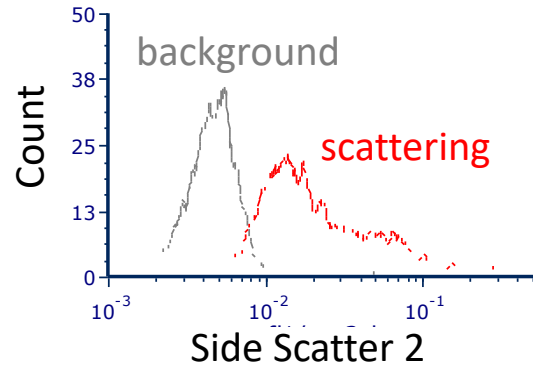
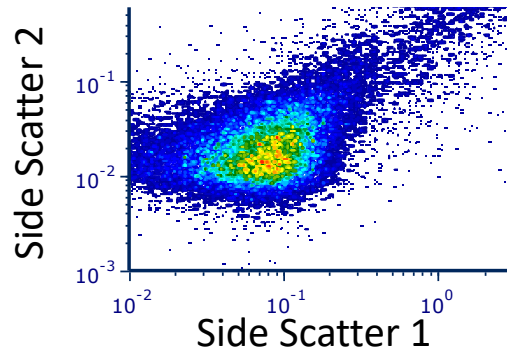


FL3 (CellBrite 650 / membrane)



- EVs: HansaBioMed HEK293 Fluo-EVs
- ZetaView: ~ 90 nm
- run A: Biotium CellBrite Steady 650, EGFP/CD63
- run B: PE/CD81

Delaware: Cell Assay: Stained VeriCell Leukocytes


lymphocytes

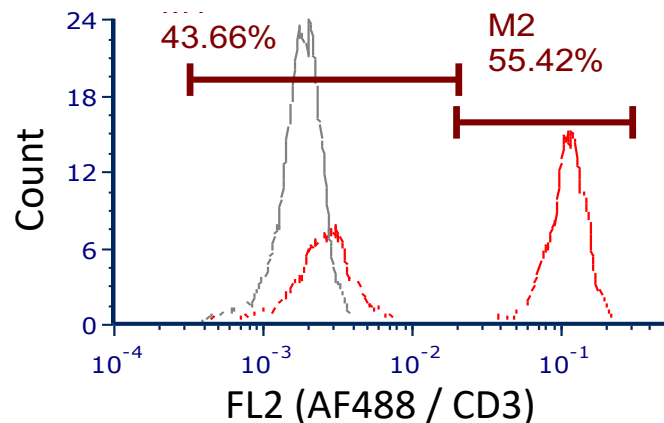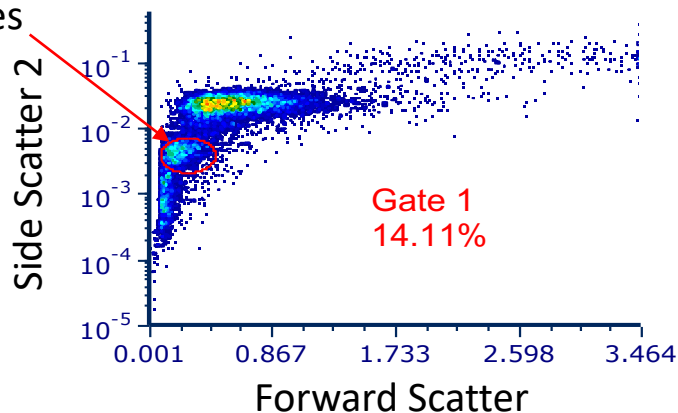

no
compensation

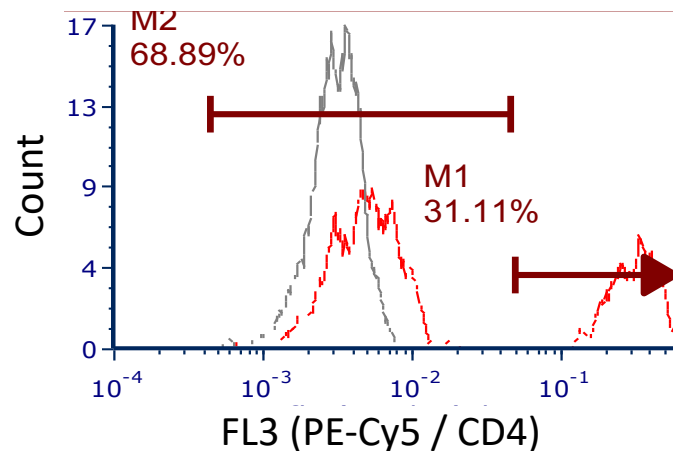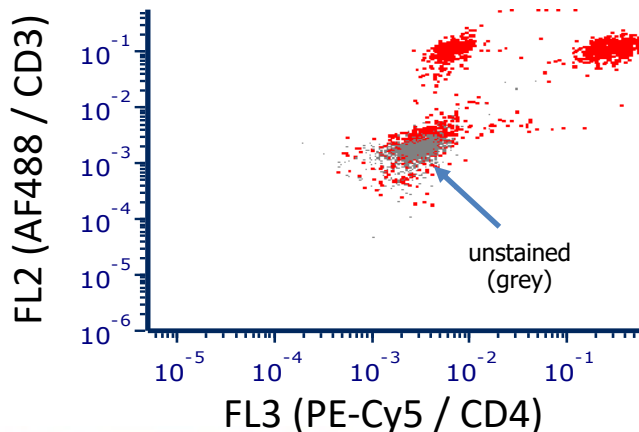

Panama Software: Instrument Control, Data Display

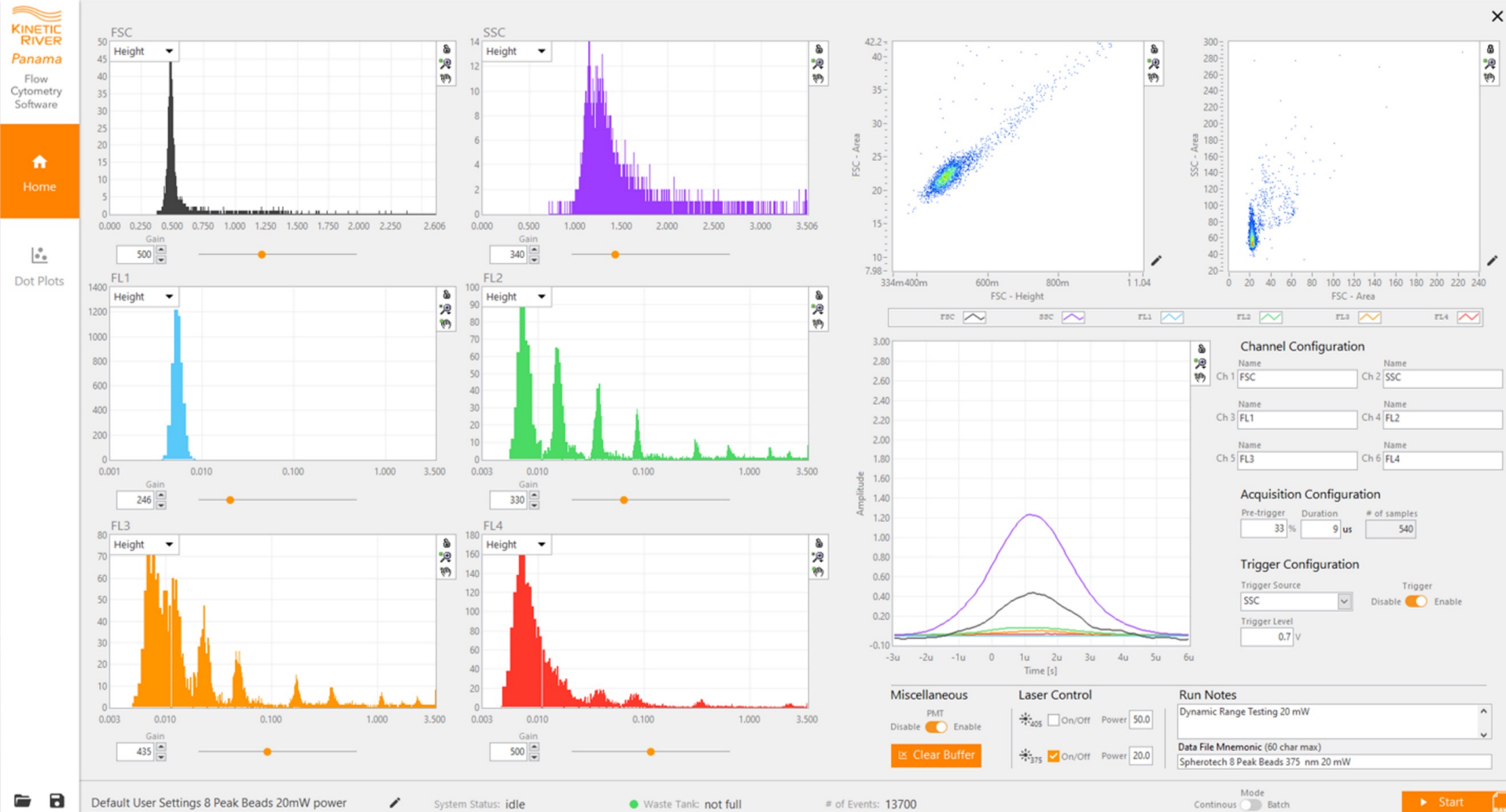

Delaware Technical Features and User Benefits

<i>Features</i>	<i>Benefits</i>
• highly stable, pressure-driven <i>Shasta</i> fluidics	→ more precise measurements
• proprietary probe wash design	→ lower background, carryover
• high-power lasers	→ higher signal-to-noise ratios
• UV/violet lasers	→ higher scattering signals
• proprietary collinear excitation architecture	→ improved scattering sensitivity
• three scattering detection channels	→ improved scattering resolution
• proprietary seven-element objective	→ improved light collection
• proprietary stray light rejection architecture	→ reduced background levels, more robust operation
• all-PMT detection	→ high sensitivity, low dark current
• proprietary signal processing algorithms	→ reduced background events
• intuitive, flexible user interface	→ easy to swap b/w different operating modes



Delaware Configurations

Basic Configuration	High Sensitivity Configuration	Five-Laser Configuration
<p>2 Lasers</p> <p>405 nm, 250 mW 488 nm, 200 mW</p>	<p>3 Lasers</p> <p>375 nm, 50 mW 405 nm, 250 mW 488 nm, 200 mW</p>	<p>5 Lasers</p> <p>375 nm, 50 mW 405 nm, 250 mW 488 nm, 200 mW 561 nm, 50 mW 640 nm, 150 mW</p>
<p>Standard Scattering</p> <p>FSC, SSC (405 and 488 nm)</p>	<p>Ultrasensitive Scattering</p> <p>FSC, SSC (375, 405, 488 nm)</p>	<p>Ultrasensitive Scattering</p> <p>FSC, SSC (375, 405, 488 nm)</p>
<p>2 Fluorescence Channels</p> <p>525/50 580/23</p>	<p>4 Fluorescence Channels</p> <p>525/50 580/23 615/24 697/58</p>	<p>6 Fluorescence Channels</p> <p>440/40 (optional) 525/50 580/23 615/24 697/58 755/35</p>

Delaware Flow NanoCytometer® for EV Analysis

- 3 configurations for broad range of needs
 - up to 5 lasers / 6 fluorescence channels
- 68-nm single-liposome detection
- 60-nm single-NP detection (PS)
- 28-nm single-NP detection (Au)
- 6-nm resolution
- 2% CVs
- HEK293, HeLa, PC3 EVs
- FL: tetraspanins, membrane, EGFP
- low background
- wide dynamic range
- up to 1,000+ events/sec
- intuitive user interface
- *also* measures cells



KineticRiver.com/Delaware
info@KineticRiver.com



KU

THE KIDNEY INSTITUTE

The University of Kansas

Toward Flow-Based Rapid Diagnostics of Kidney Disease.

Christopher J. Ward, MBChB, PhD

Division of Nephrology and Hypertension

Jared Grantham Kidney Institute

University of Kansas Medical Center

030724

Speaker Bio

- Trained as a medical doctor 1981-86 (Edinburgh, Scotland).
- PhD Birmingham University UK 1986-1990.
- Worked on ADPKD from 1991-present (33 years).
- Involved in the cloning of *PKD1*, *TSC2* (Oxford) and *PKHD1* (Mayo Rochester).
- Reagent development (monoclonal antibodies).
- Exosome hypothesis of ADPKD.
- Development of assays for diagnosis and prognosis of ADPKD.


Autosomal dominant polycystic kidney disease (ADPKD)



About PKD

- Affects 1:800 people, 425,000 in the US.
- Dominant disease.
- Mainly private mutations.
- Long therapeutic window.
- Multiple treatment in the pipeline, including strategies to increase the gene expression.
- Need for a rapid, low cost and non invasive monitoring test, better than MRI.

Autosomal dominant polycystic kidney disease (ADPKD)

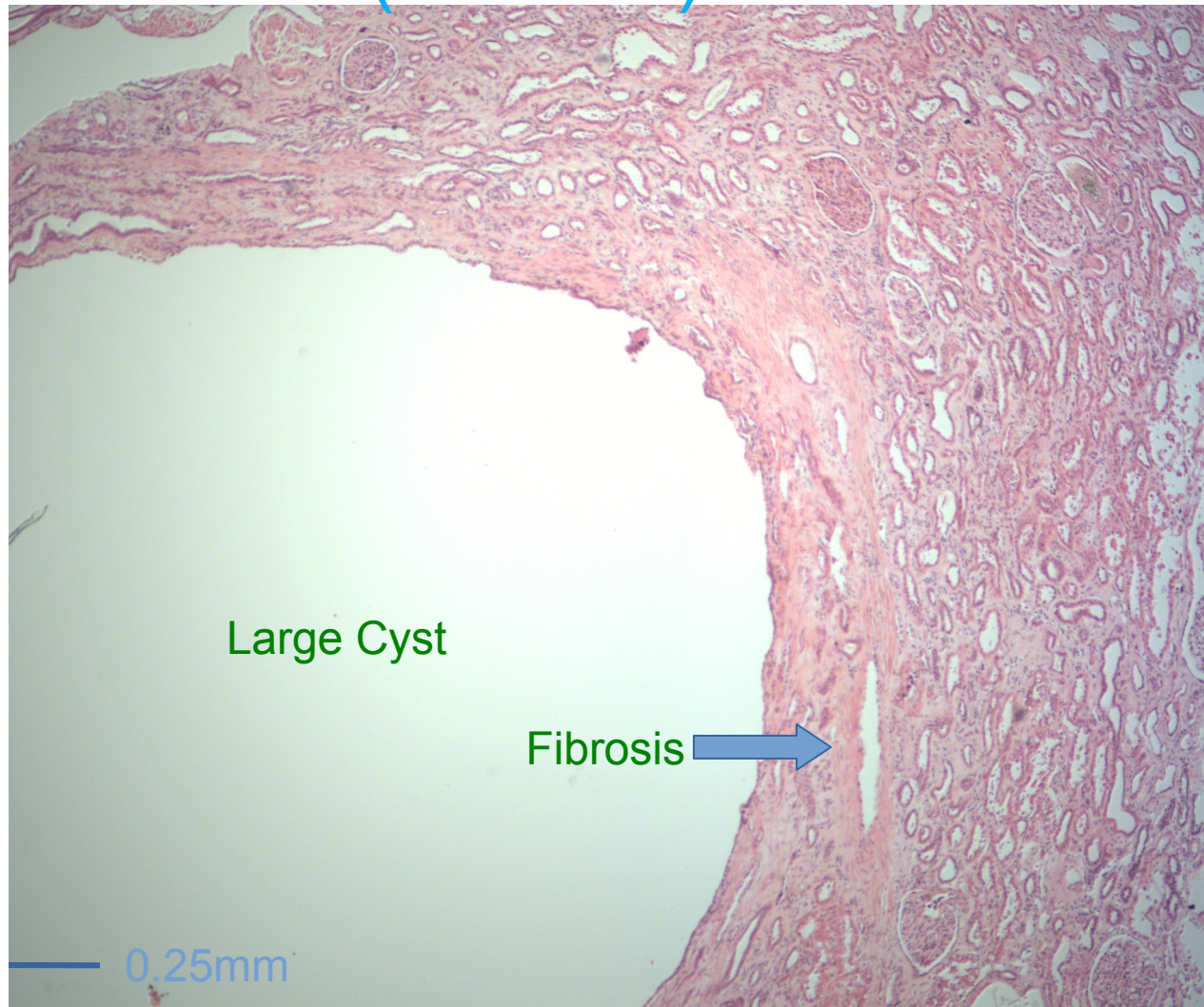


Kidney from normal (WT) and ADPKD mice

RED:

S-phase


GREEN:

Apoptotic

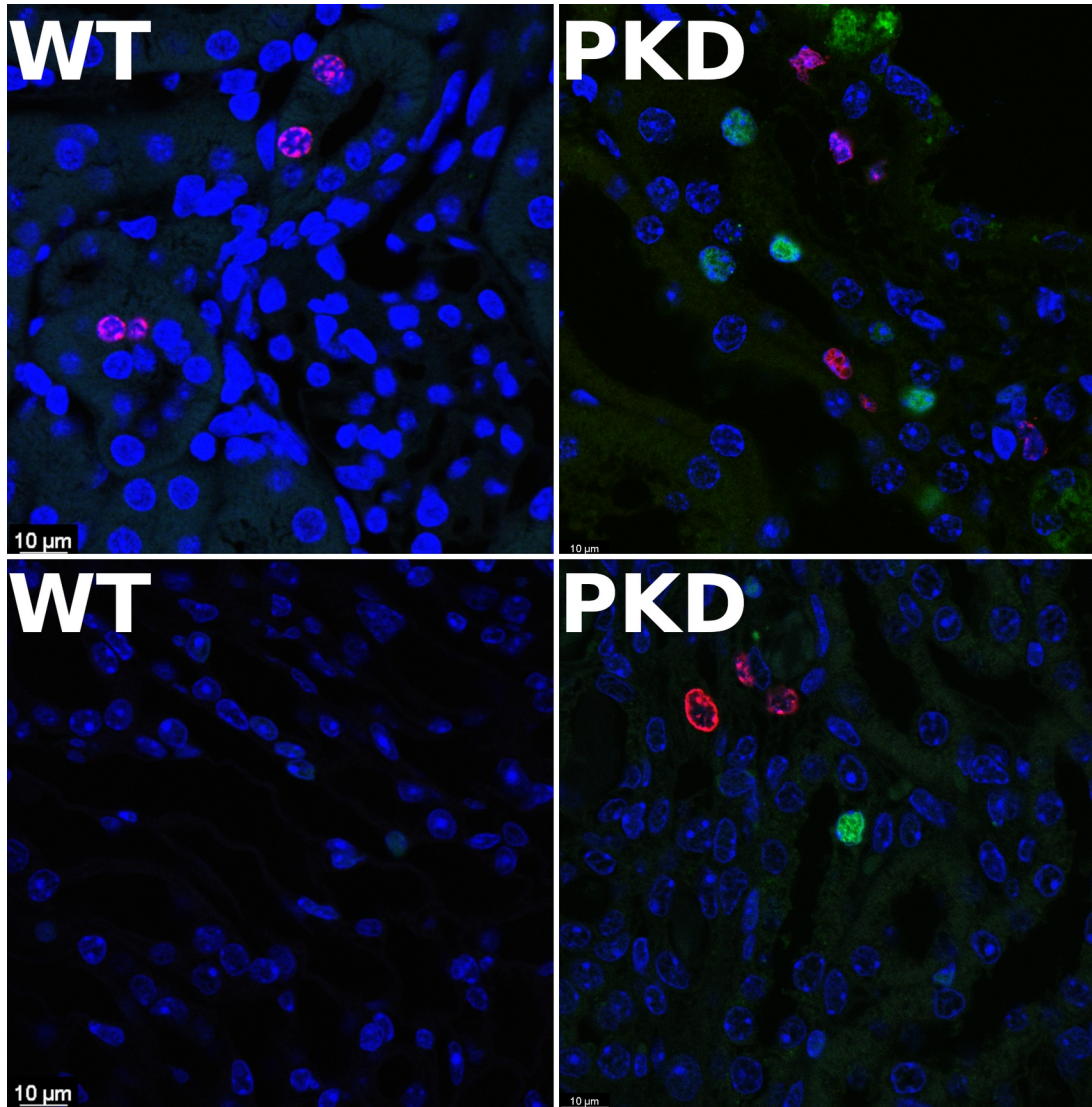


Polycystic kidney disease genes

We were responsible for identifying the PKD1 gene (1994)

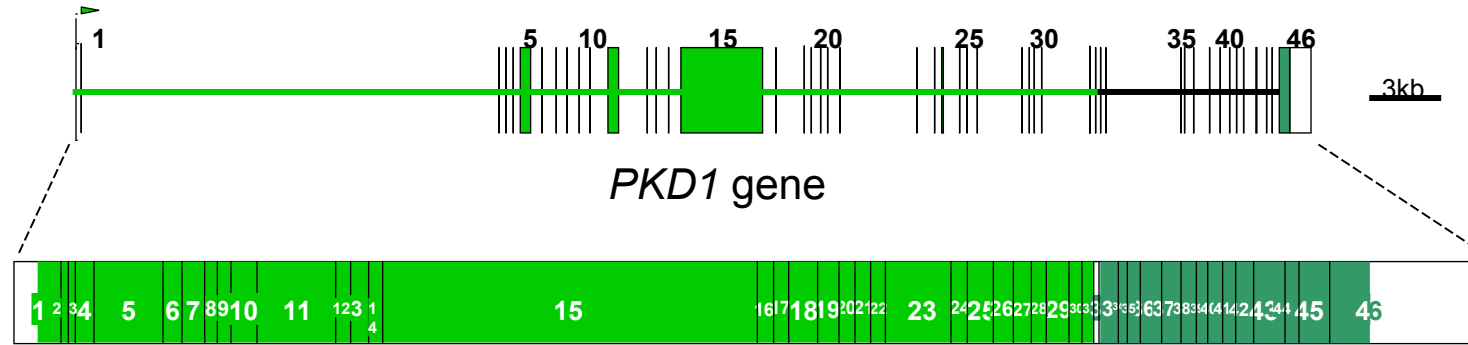

PKD1 ~85% cases 16p13.3

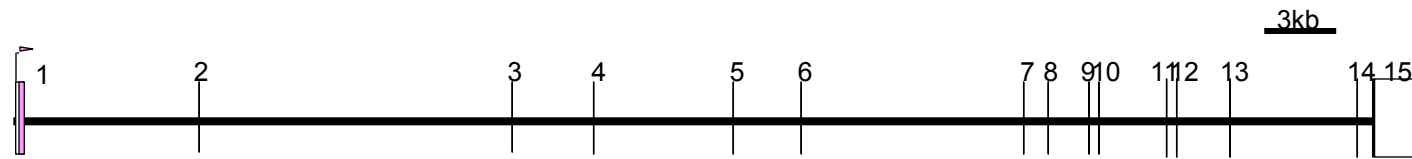


PKD2 ~15% cases

4q21-q23


European PKD1 Consortium, Ward CJ *et al.* Cell 1994

We also developed a model of the proteins spanning the membrane and developed monoclonal antibodies to the proteins.

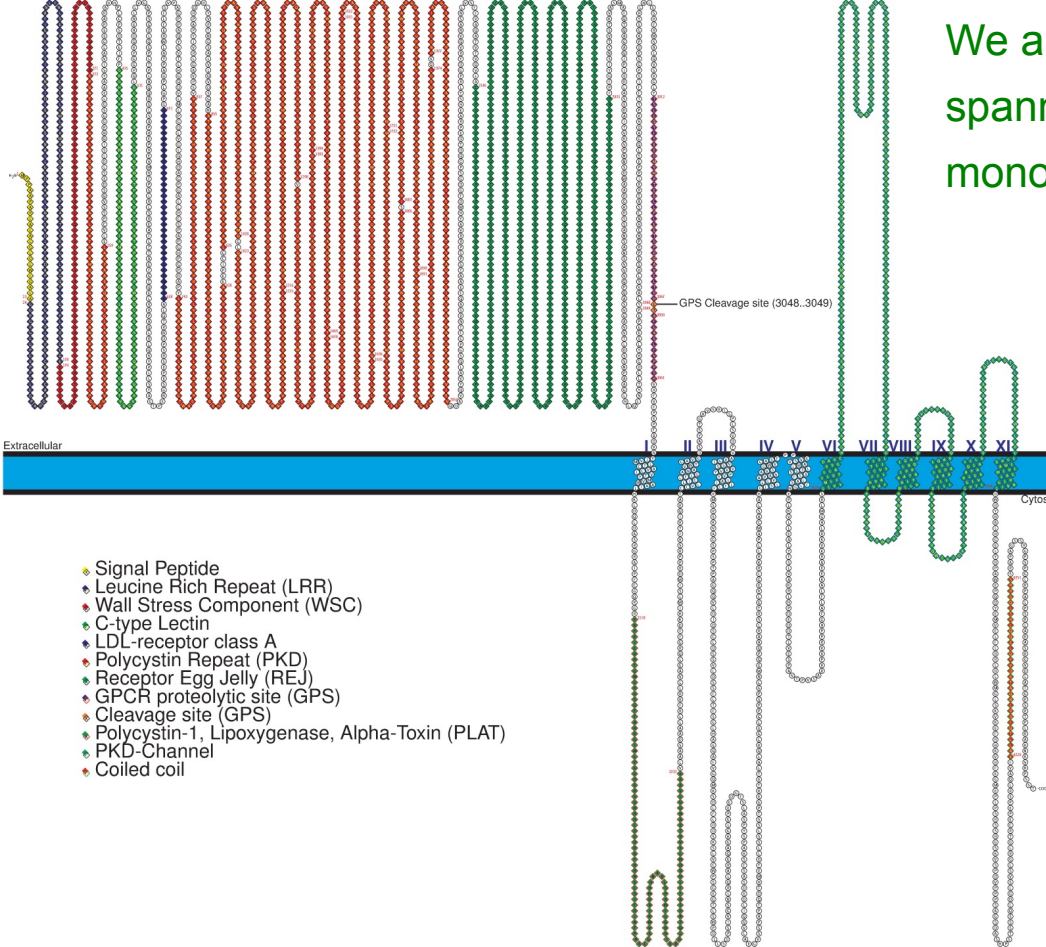


- ◆ Signal Peptide
- ◆ Leucine Rich Repeat (LRR)
- ◆ Wall Stress Component (WSC)
- ◆ C-type Lectin
- ◆ LDL-receptor class A
- ◆ Polycystin Repeat (PKD)
- ◆ Receptor Egg Jelly (REJ)
- ◆ GPCR proteolytic site (GPS)
- ◆ Cleavage site (GPS)
- ◆ Polycystin-1, Lipoxigenase, Alpha-Toxin (PLAT)
- ◆ PKD-Channel
- ◆ Coiled coil

Polycystin-1
4300aa

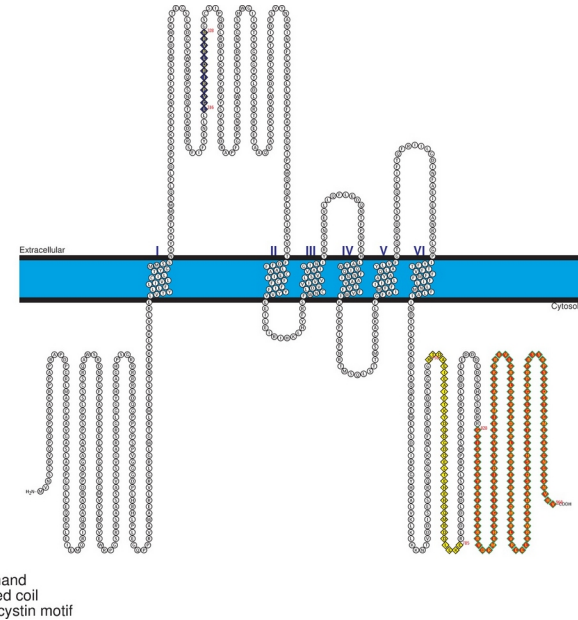

Polycystin-2
1000aa

PC1 enriched urinary exosomes are 100-150nm diameter

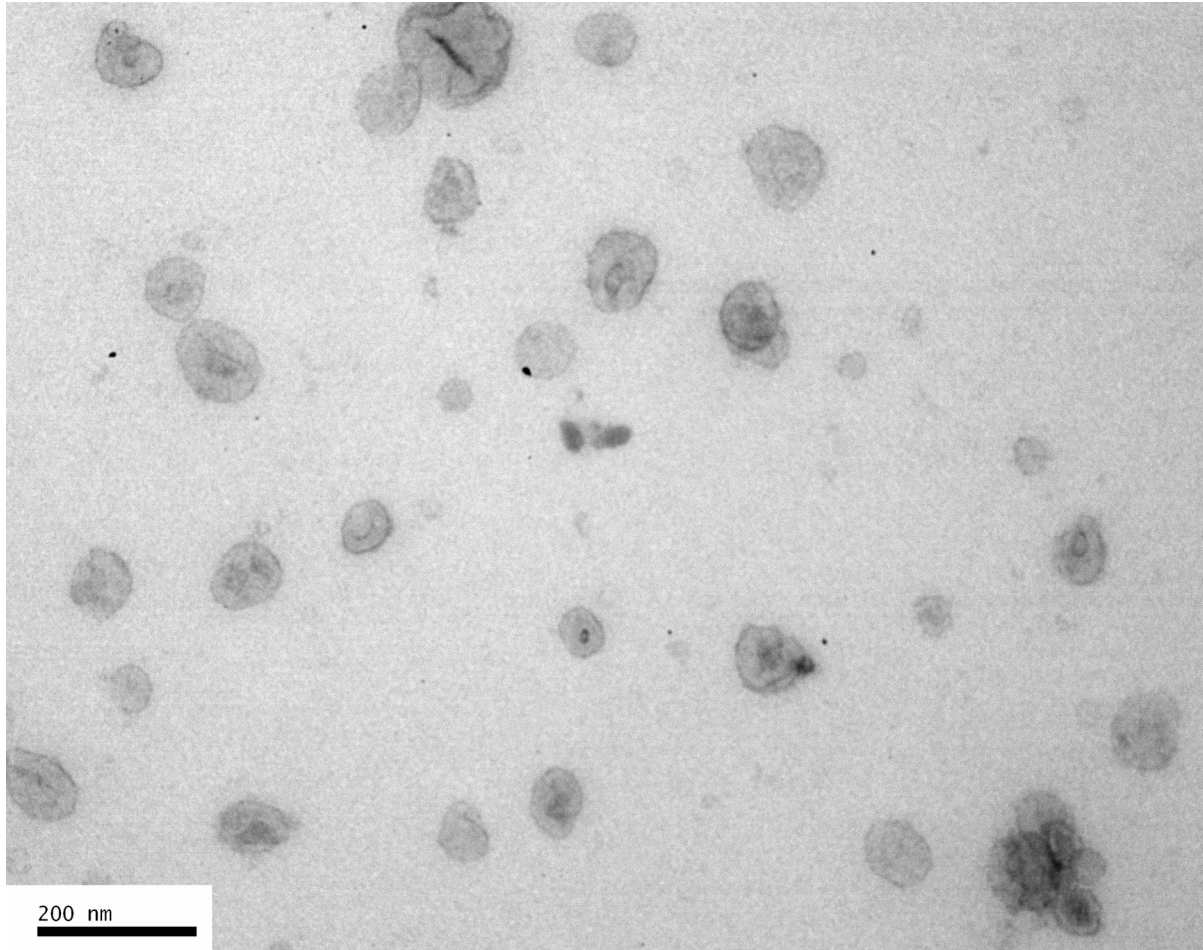


Purified from human urine


Volcano Plot of *PKD1* vs Normal

Increase statistical significance

$-\log_{10}$ (adjusted p-value) against \log_2 (fold change PKD1 vs WT)

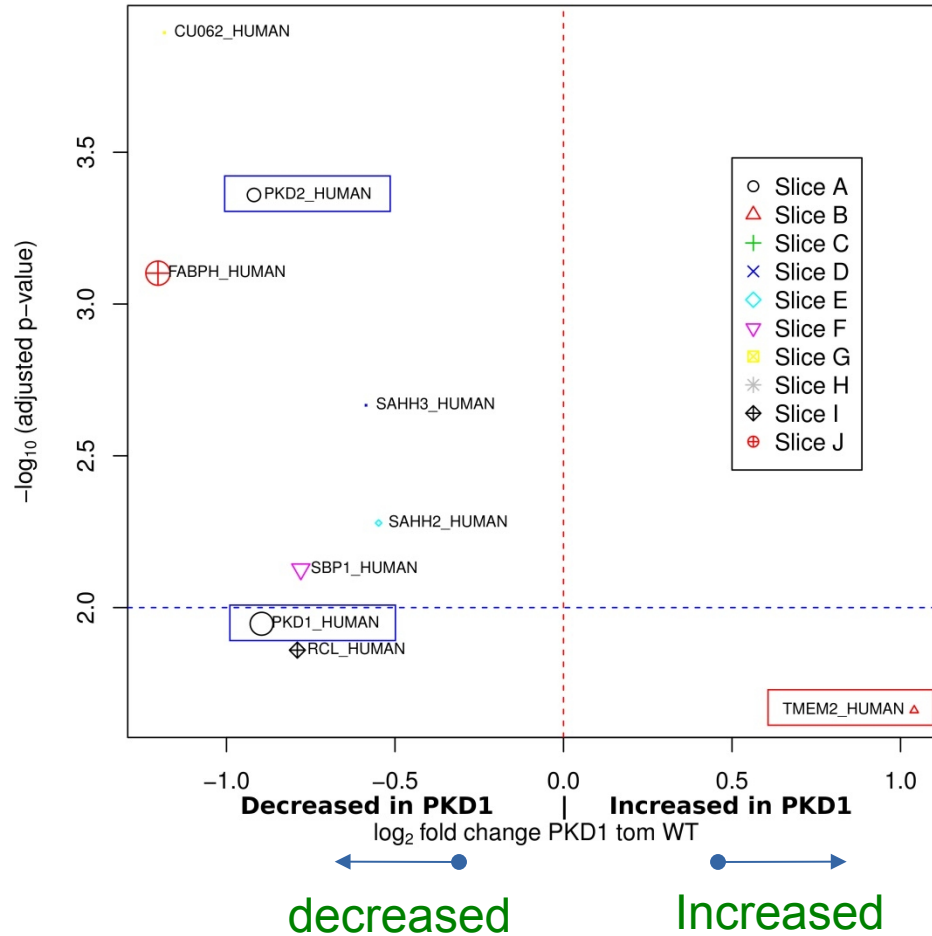

Western blot analysis of PC1 levels controlled by CD133.

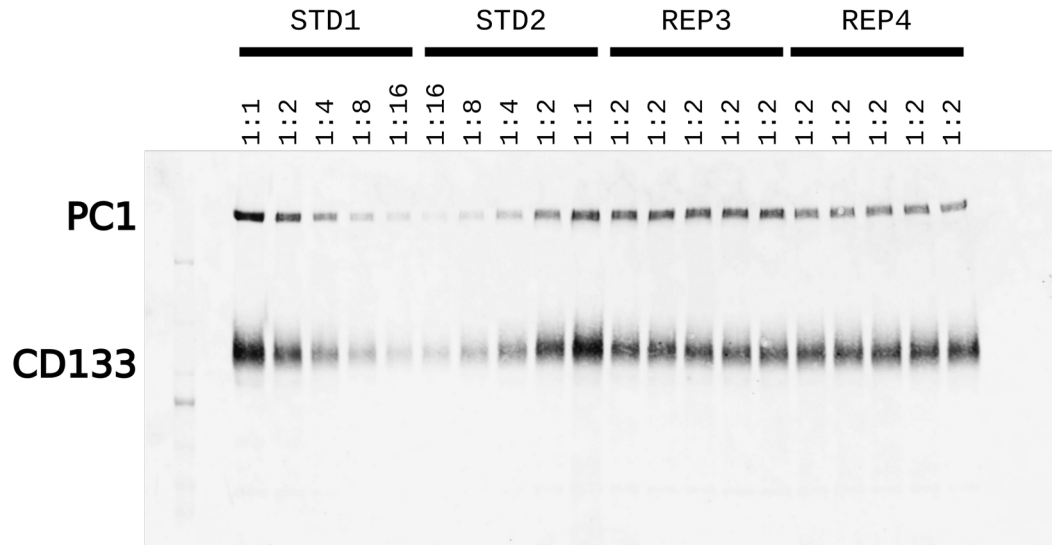


Linear dose response in batch CV 12%

This Western shows a titration of EVs STD1 & 2 and two repeat loadings to determine the variation in the assay.


A Western blot can be used to diagnose ADPKD but the variability inherent in the assay is such that we cannot determine severity or prognosis.

PC1/CD133 ratio is a promising biomarker for ADPKD.

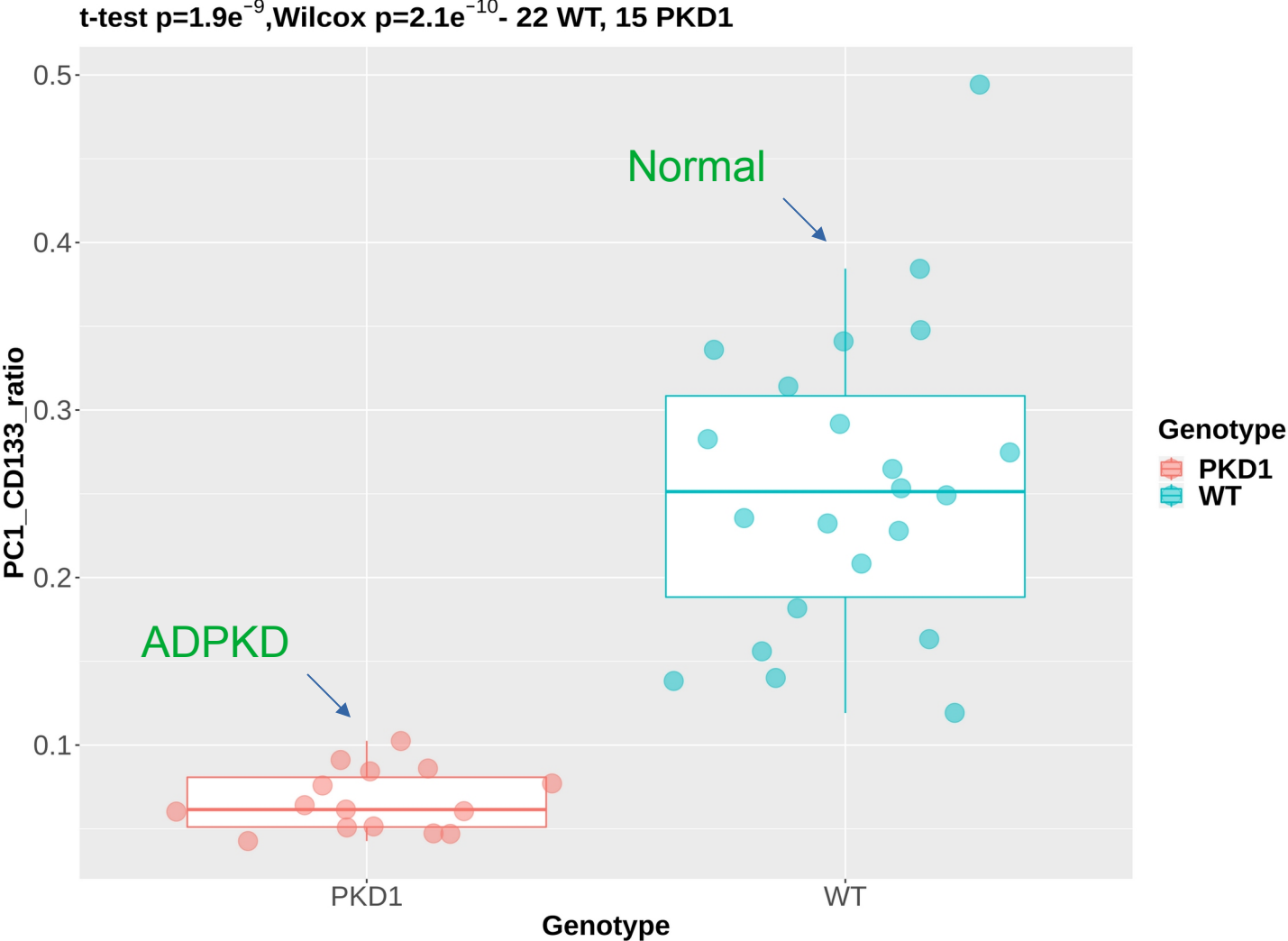


MS/MS data HtTKV against PC1_CD133 ratio, $R^2=0.63$

This shows that urine EVs contain important information about disease severity.


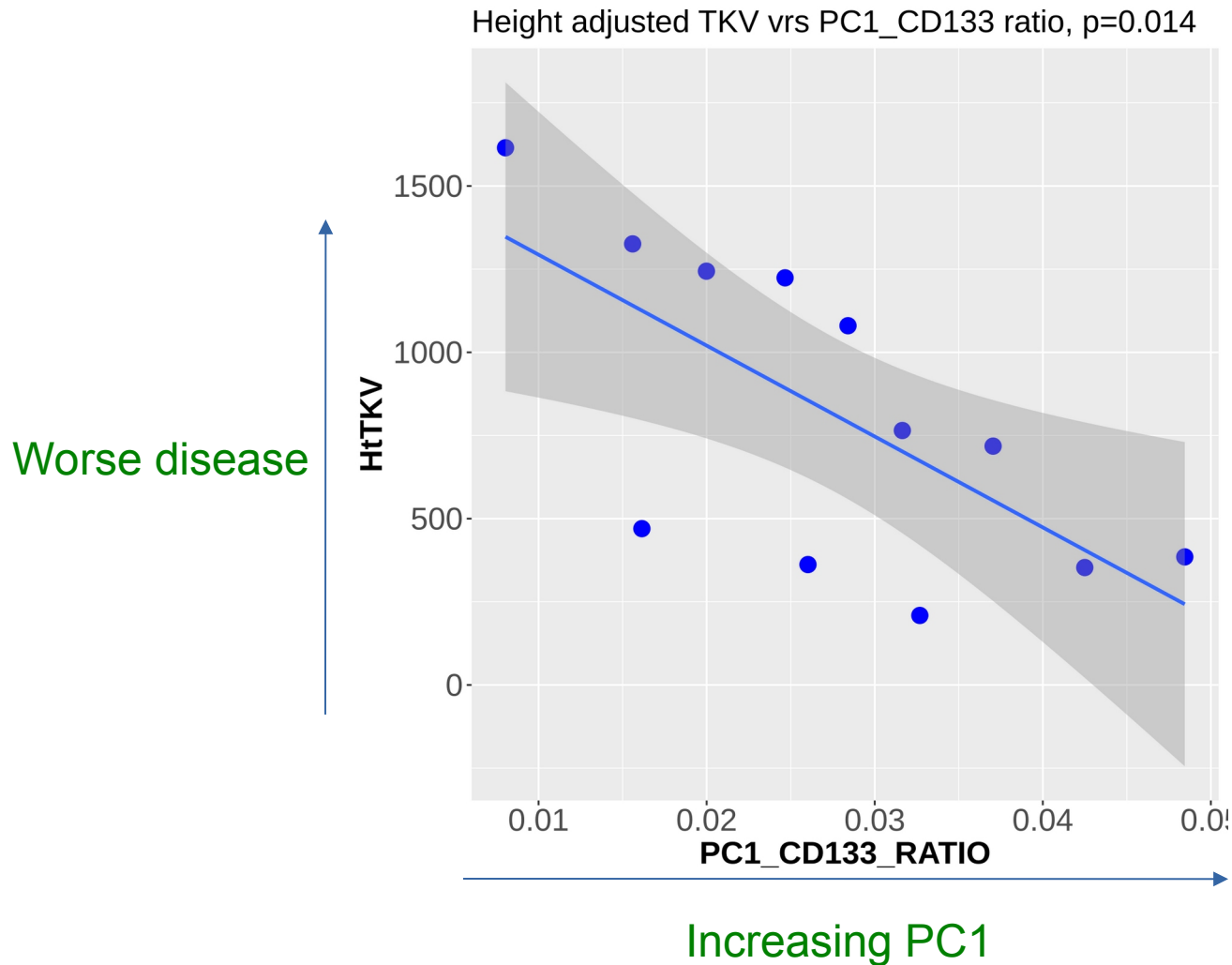


The anti-PC1 monoclonal antibody can detect urinary exosomes.


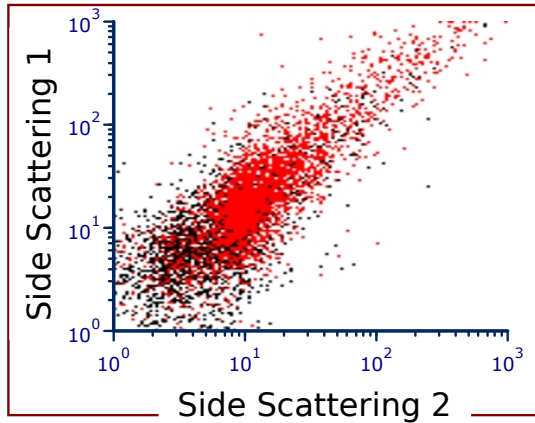


Results collected on the *Delaware* Flow NanoCytometer

no membrane stain

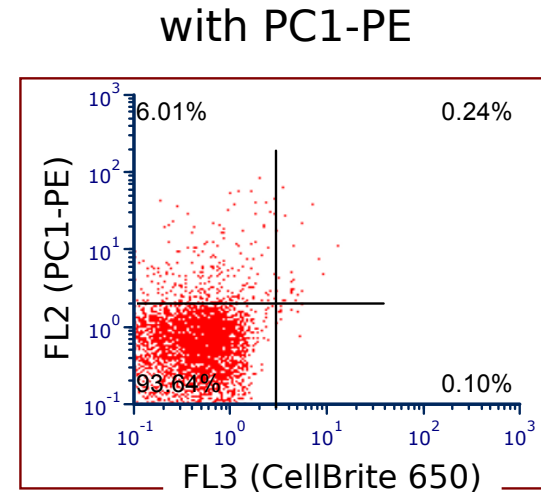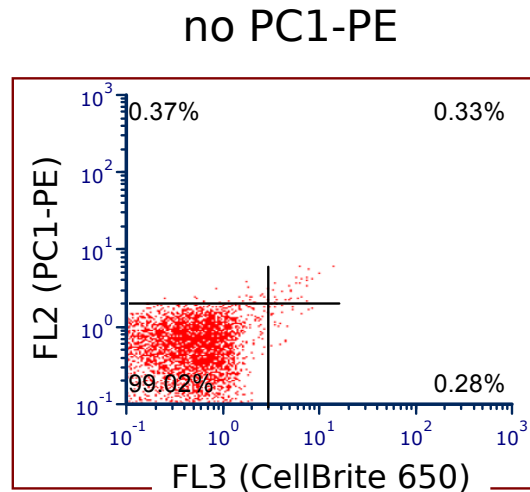

with membrane stain

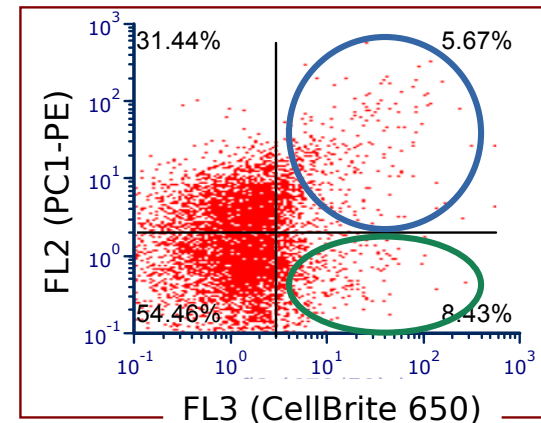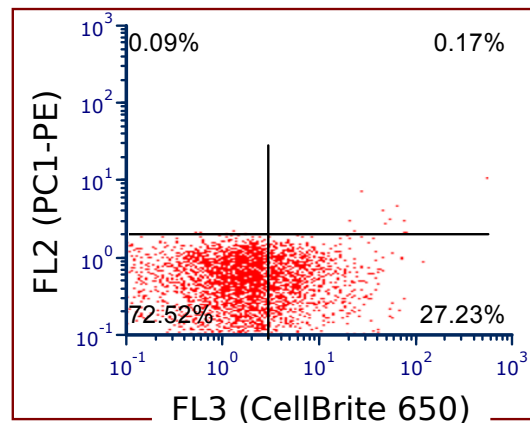

Other possibilities:

Proteins decreased in ADPKD, for which we have mAbs:

- Polycystin-2 (PC2)
- Exosomal polycystin-1 interacting protein (EPCIP)
- *C16orf89* protein
- Fibrocystin

There is one protein that is increased in ADPKD:

- TMEM2 (CEMIP2) work in progress.

Acknowledgements

- *Christopher J. Ward*
- *Wendy Lea*
- *Lesya Zelenchuk*
- *Madhulika Sharma*
- *Kerri McGreal*
- *Darren Wallace*
- *Gail Reif*

